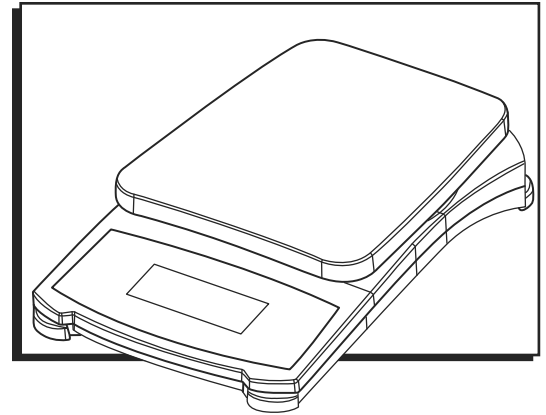
ULINE H-8109, H-8110
 H-8111

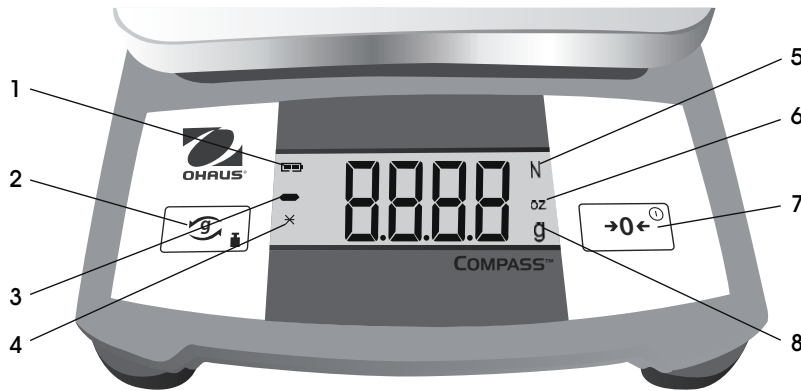
1-800-295-5510
 uline.com

OHAUS CX COMPASS SCALE

OVERVIEW OF CONTROLS

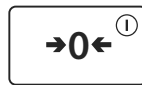
CONTROL PANEL AND DISPLAY

CONTROL PANEL PARTS

#	DESCRIPTION
1	Battery Indicator
2	Unit Cal Menu Button
3	Negative Sign
4	Stability Indicator
5	Symbol for Newtons
6	Symbol for Ounces
7	On/Off Zero Button
8	Symbol for Grams

CONTROL FUNCTIONS

BUTTON		
PRIMARY FUNCTION (Short Press)	Changes unit of measure.	Turns the scale on. If the scale is on, sets zero.
SECONDARY FUNCTION (Long Press)	Enter user menu.	Turns off the scale.

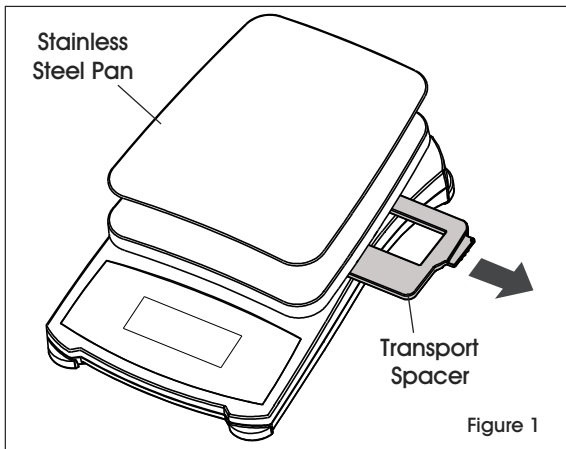
SETUP

WARNING! Verify that the AC adapter input voltage matches the local AC power supply. Use the scale only in dry locations. For best results, use the scale in the specified optimum operating temperature. Do not operate the scale in hostile or unstable environments. Use the scale with care, gently loading items to be weighed onto the center of the pan. Avoid rough treatment to prolong life of scale. Indoor use only.

REMOVING TRANSPORT SPACER (H-8109 ONLY)

1. To remove the red transport spacer, pull it out from the side of the scale. (See Figure 1)

INSTALLING THE PLATFORM

1. Place the stainless steel pan onto the weighing platform. (See Figure 1)

CONNECTING POWER

BATTERY

1. Remove the battery cover on the bottom of the scale.
2. Place three AA batteries into the compartment.

CAUTION! Do not use excessive force or press on the weighing pan.

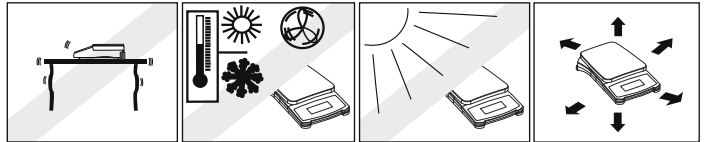
3. Re-insert the battery cover.

AC POWER (OPTIONAL. H-8112 SOLD SEPARATELY.)

1. Connect the AC adapter to the receptacle at the rear of the scale.
2. Plug the AC adapter into a properly grounded electrical outlet.

SELECTING THE LOCATION

1. Place on a firm, steady surface with sufficient space.
2. Avoid excessive air currents, vibrations, heat sources or rapid temperature changes.

OPERATION

POWER UP TEST

When the scale is turned on, all display segments will be displayed for a few seconds, indicating that the unit is self-adjusting to zero. Once 0 is displayed, the scale is ready for use.

STABLE READING INDICATION

A star indicator, *, will appear in the lower left corner of the display when a stable reading has been reached.

OVERLOAD

If the applied load exceeds the capacity of the scale, E will appear on the display and the load should be removed immediately. The scale will return to normal operation.

TARE FUNCTION

Values can progressively be added to a sample. By pressing the **ON/OFF ZERO** button, the scale display returns to zero. After adding additional mass, press **ON/OFF ZERO** to zero the scale again. Additional mass may be added up to the capacity of the scale.

NEGATIVE VALUE

When a load is removed from the scale, any zeroed value will be displayed as a negative number. To return to normal operation, the zeroed value can be cancelled by pressing **ON/OFF ZERO** button.

OPERATION CONTINUED

LOW BATTERY INDICATION

The display will show *L0* when the batteries are weak and need to be replaced.

AUTO SHUT-OFF

To extend battery life, the scale will automatically turn off after approximately four minutes if no active weighing is occurring. This feature is active with battery operation only.


CALIBRATION

For best results, calibrate the scale at regular intervals, especially if using for prolonged periods. Calibration weights are not provided with the scale.

CALIBRATION VALUES

MODEL #	CALIBRATION POINT (g.)
H-8109	200
H-8110	1,000
H-8111	2,000

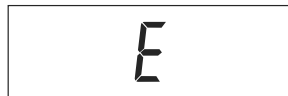The message *E* will appear if the calibration steps are not followed or the wrong weight was used.



1. Press and hold the button until *MENU* is displayed. Release the button, first menu *CAL* will be displayed.

NOTE: The calibration process can be aborted by pressing .

2. Press to start span calibration.

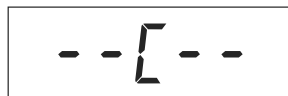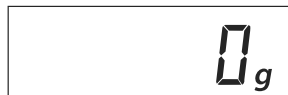3. The display shows *0*. With no weight on the pan, press the button to establish the zero point. The display shows *--[--* while the scale stores the zero load signal.


4. The display shows the span calibration point, for example, 200g. Place the specified calibration weight on the pan and press the button. The display shows *--[--* while the scale stores the calibration point signal.

5. The display shows *[-done-]* and returns to the normal weighing mode.

BACKLIGHT

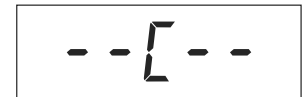The backlight settings can be adjusted using the menu. Press and hold the button until *MENU* is displayed. Release the button, first menu *CAL* will be displayed. Press to select desired option.

MENU OPTION	SETTING
<i>b.Auto</i>	Scale turns on light automatically with any button or weighing operation.
<i>b.ON</i>	Backlight is always on.
<i>b.OFF</i>	Backlight is always off.

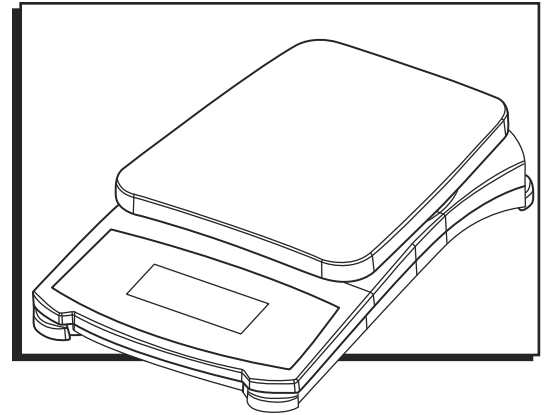
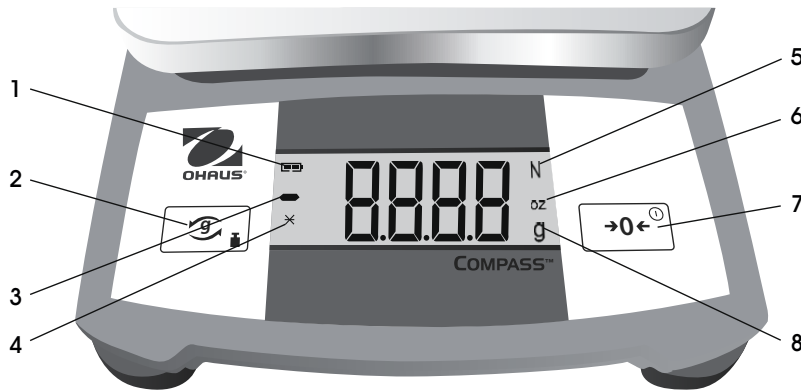
Once desired setting is displayed, press to select and return to current weighing mode.

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